

## PERMAGRAPH® C

for the computer controlled  
measurement of magnetization curves  
of hard magnetic materials



### • Introduction

The PERMAGRAPH® is the world-wide known standard equipment for the measurement of magnetic characteristics of permanent magnets in research, development, quality control and production.

The PERMAGRAPH® C is an automatic computer-controlled precision version of this measuring instrument. Equipped with a dual-channel Electronic Fluxmeter EF 7 and the high-performance software package PERMA, it facilitates quick and reliable measurements. The PERMAGRAPH® C will meet your present and future requirements. Our instruments are known for their quality and long lifetime.

The system design conforms to the standards IEC 60404-5, ASTM A977/A977M and DIN IEC 60404-5 (replacement for DIN 50470, DIN EN 10332). The temperature upgrade conforms to IEC 61807TR and DIN IEC 68/190/CDV. The equipment has been optimized over many years and has been made specifically to meet the requirements of our customers.

Among others, the following measurements are possible:

- Automatic measurement of the hysteresis loops of permanent magnets
- Determination of characteristic quantities like remanence, coercivity, maximum energy product, etc.
- Measurements at temperatures up to 200 °C with heating poles and temperature surrounding coils
- Measurements on the surface of magnets using pole coils

MAGNET-PHYSIK has 75 years of experience in making measuring instruments for the characterization of magnetic materials.

## • Applications



Depending on the model of the PERMAGRAPH® C and its particular accessories the following measurements are possible:

**Measurements on AlNiCo, Ferrite or Rare Earth magnets (e.g. Sm-Co or Nd-Fe-B)**  
using J-compensated surrounding coils.

**Important:**

For the saturation of rare earth magnets an impulse magnetizer and a magnetization coil are required.

**Measurement of AlNiCo, Ferrite or Rare earth magnets at high temperatures**

with up to 200 °C heatable poles and temperature resistant J-compensated surrounding coils.

You can find technical explanations and further topics, e.g. how hard magnetic powders are measured with the PERMAGRAPH®, in our booklet

**MAGNETIC MEASURING TECHNIQUES**

by Dr. Erich Steingroever  
and  
Dr. Gunnar Ross

It can be obtained free of charge upon request (also available as a pdf file).

**Measurements on Ferrite magnets**


using measuring poles with embedded pole coils.

**Measurement of Ferrite segment magnets**

with segment poles with incorporated pole coils, the poles are shaped corresponding to the radii of the segment magnet.

**Magnet shapes**

Examples of specimen shapes that can be measured with pole coils in the PERMAGRAPH® C without the need of determining their cross section:

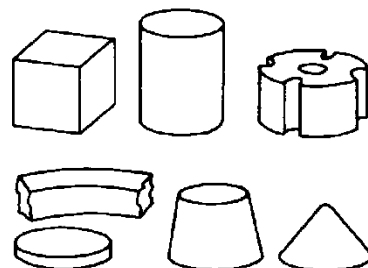

## • Product Family



The PERMAGRAPH® C can be composed from a variety of components and accessories.

For the most common measurement tasks we offer packages which can be completed by further components according to your application.

On the following pages the components that are included in the packages are listed. You will also find descriptions and technical details of all available accessories.

All components are calibrated in our qualified calibration laboratory. The calibration is traceable to national standards. A proprietary calibration certificate is included in the delivery.

## • Specialties of the MAGNET-PHYSIK Measuring Method

For measuring the field strength  $H$  we use either integrated or separate field strength measuring coils. Therefore, a Hall effect gaussmeter is superfluous. When Hall sensors are used to measure  $H$  next to the specimen, additional space for the probe and an additional gaussmeter are required. Hall sensors are very fragile and can easily be damaged. Because of the linearity error and the temperature dependence of the sensitivity of a Hall sensor, corrections are necessary to achieve an acceptable accuracy. Additional errors can arise from the facts that the Hall probe has always to be aligned truly perpendicular to the magnetic field direction and that it is, due to the small active area, more sensitive to local field strength variations. Therefore, it is the better solution to use a field strength measuring coil instead of a Hall probe.

If a surrounding coil is used, the field strength and the polarization are measured with only one combined coil system. As no space for the Hall probe is required, specimens of a thickness down to only 1 mm can be measured.

• Packages, Upgrades and Options

<p><b>PERMAGRAPH® C - 400</b> for measurement at ambient temperature</p> <ul style="list-style-type: none"> <li>1 Electronic fluxmeter EF 7 with two integrator modules FM 1</li> <li>1 Power supply SVP 5</li> <li>1 Cabinet for Permagraph® C - 400</li> <li>1 Electromagnet EP 3</li> <li>2 Poles P 0/0 - ø 92 mm</li> <li>1 J-compensated surrounding coil JH 26-1 (other diameter on request)</li> <li>1 Computer</li> <li>1 Software PERMA for Windows</li> <li>1 Training in our works</li> </ul>	<p><b>PERMAGRAPH® C - 400</b> for measurements at temperatures up to 200 °C</p> <ul style="list-style-type: none"> <li>1 Electronic fluxmeter EF 7 with two integrator modules FM 1 and one temperature module TM 1</li> <li>1 Power supply SVP 5</li> <li>1 Cabinet for Permagraph® C - 400</li> <li>1 Electromagnet EP 3</li> <li>2 Poles P 0/0 - ø 92 mm</li> <li>1 J-compensated surrounding coil JH 26-1 (other diameter on request)</li> <li>2 Heating poles PT 200</li> <li>2 Thermocouple</li> <li>1 Temperature surrounding coil JHT 40-4 (other diameter on request)</li> <li>1 Computer</li> <li>1 Software PERMA for Windows</li> <li>1 Training in our works</li> </ul>
<p>A PERMAGRAPH® C - 400 can also be upgraded later for measurements at temperatures up to 200 °C.</p> <p>The temperature module TM 1 is computer-controlled. The desired temperature is set from the computer and the measurement temperature is automatically saved with the measurement.</p>	<p><b>Upgrade package TM 1</b> for PERMAGRAPH® C - 400 for measurements at temperatures up to 200 °C</p> <ul style="list-style-type: none"> <li>1 Precision temperature control unit TM 1 with independent control for the two heating poles. The TM 1 is installed into the EF 7.</li> <li>2 Heating poles PT 200</li> <li>2 Thermocouples</li> <li>1 Temperature surrounding coil JHT 40-4 (other diameter on request)</li> </ul>
<p><b>Option Electromagnet EP 5</b> EP 5 to be supplied <i>instead of</i> the electromagnet EP 3 which is included in the packages above</p>	
<p><b>Option PT 200 FeCo</b> Heating poles PT 200 FeCo to be supplied <i>instead of</i> the PT 200 which are included in the upgrade packages</p>	
<p><b>Detailed information on the individual components is given on the following pages.</b> <b>For further accessories see page 11 and following.</b></p>	

## Technical Data

<b>Cabinet for PERMAGRAPH® C - 400</b> Houses a controller for the electromagnet power supply and polarity reversal, and the electrical distribution.	Width	:	515 mm
	Height	:	300 mm
	Depth	:	620 mm
	Weight	:	20 kg
	Mains standard*	:	3 x 380 - 415 V
		:	50 - 60 Hz
	Mains optional	:	1 x 230 V
		:	50 - 60 Hz

<b>Electronic Fluxmeter</b> To measure the polarization $J$ as well as the field strength $H$ (Dual-channel version)	<b>EF 7</b>	Detailed technical data and a description of the properties of this instrument can be found on one of the next pages. Weight:	:	3.2 kg
<b>PERMAGRAPH® Power Supply</b> Power supply for the electromagnet, with minimum ripple voltage, electronic remote control	<b>SVP 5</b>	Output voltage Output current Output power Mains*  Mains optional  Weight	:	0 - 120 V 0 - 25 A 3 kW 3 x 380 - 415 V 50 - 60 Hz  1 x 230 V 50 - 60 Hz  14 kg
<b>PERMAGRAPH® Electromagnet</b> For the magnetic field excitation to record magnetization curves of hard magnetic materials. The pole distance is manually adjustable.	<b>EP 3</b>	Max. field strength 92 mm standard poles, 2 mm air gap Bottom pole size Standard pole diameter Air gap (without poles) Air gap (with standard poles) Nominal short-time electrical load Depth x width x height Weight	:	approx. 1700 kA/m, (21.5 kOe) 100 mm x 100 mm 92 mm 73 - 102 mm 0 - 63 mm 3 kW (310 x 500 x 600) mm <sup>3</sup> 128 kg
<b>PERMAGRAPH® Electromagnet</b> For the magnetic field excitation to record magnetization curves of hard magnetic materials. The pole distance is manually adjustable.	<b>EP 5</b>	Max. field strength 92 mm standard poles, 2 mm air gap 65 mm FeCo poles, 2 mm air gap Max. pole diameter Air gap, without poles Air gap, with standard poles Nominal short-time electrical load Depth x width x height Weight	:	approx. 2200 kA/m, (27.5 kOe) approx. 2500 kA/m, (32 kOe) 92 mm 0 - 130 mm 0 - 110 mm 3 kW (430 x 480 x 830) mm <sup>3</sup> 220 kg

\*Other line voltage configurations may require an internal or external transformer. Please contact us.

## • Operation



The PERMAGRAPH® can automatically record the demagnetization curves of permanent magnets. The measuring method is standardized and verified by round robin tests.

Central component of the measuring instrument is an electromagnet that is used to magnetize and demagnetize the magnet specimen in order to cycle through the hysteresis loop. The magnetic field strength  $H$  and the magnetic polarization  $J$  are measured simultaneously with special measuring coils. Two integrators (fluxmeter channels) are required to process the output signals of the coils. The measuring coils and their parameters are automatically identified by the fluxmeters. The PERMAGRAPH® C uses an EF 7 precision fluxmeter with unique digitally compensated integrators.

The two fluxmeter channels are synchronously triggered. They are operated from a computer. The measured data are received from the fluxmeter via the USB interface and shown on the computer display in real-time. This allows checking the results already while the measuring procedure is in progress. The measuring process is controlled via a digital computer interface.

The computer also allows an evaluation and output of the measuring data, so that the measuring results are directly available.

Usually only the second quadrant of a hysteresis loop, the demagnetization curve, is of interest for the application of permanent magnets. Therefore, the measurements can be automatically stopped after the coercivity  $H_{cJ}$  is reached.

The measurement is carried out quasi-statically, which means that the demagnetization curve is recorded rather slowly, avoiding eddy current problems.

The specimen and measuring parameters can be stored. Thus, in the case of new measurements only a few new inputs must be made.

## • Features of *PERMA* – Software for *PERMAGRAPH*®

- Flexible, user friendly operation using menus, function keys and shortcut keys or icons
- Extensive help files, context-sensitive help
- Convenient input of measuring parameters
- Saving and opening of parameters and measuring data
- Multiple measurements can be simultaneously opened for easy comparison
- Existing measurements can serve as templates for new measurements
- Automatic calculation of measured results
- Automatic saving of measured data, parameters and results (e.g. under a test name or number)
- Saving of a group of associate measurements into a single file
- Export of measuring data, parameters and results to text files or to Microsoft Excel® files
- Export of parameters and results to text databases or SQL databases, a database viewer is integrated in the software
- Print preview for measurement diagrams with curves, parameters and results
- Output of measurement diagrams to a printer or pdf writer (printer and pdf writer not included)
- Copying of measuring diagrams and result lists via the Windows® clipboard
- Saving measuring diagrams as pictures (bmp, gif, jpeg, png) for easy distribution
- Various possibilities for customer specific output design as selection of curves, calculated results, units, measuring parameters, user-definable information texts, company logo, etc.
- User-definable limit classifiers for all results (out of range results are shown in red or boldface)
- Display of multiple curves in one diagram including results
- Optional display of averaged results of multiple measurements
- Different output diagram layouts can be saved for easy switching between different outputs
- Selectable units for magnetic quantities, temperatures, specimen dimensions and other parameters
- Full support of SI and CGS units in software and output, changing units is possible at any time
- Simultaneous display of results in mixed SI and CGS units can be configured by the user
- The number of significant digits to be displayed for results can be selected by the user
- Creating of result lists for multiple measurements incl. the possibility of saving, copying and printing
- Program menu access can be restricted for selected users (password protection)
- Language separately selectable for program menus and output (English, German, French, Spanish, Polish, Czech, Slovak, Russian, Chinese (simplified))
- Microsoft Windows® 10 / 11 compatible

## • Parameters

- Default parameters minimize the number of necessary inputs
- Automatic measuring coil type and coil data identification
- Calculation of the cross-sectional area from specimen geometries
- Input of measurement and specimen identification data to predefined or user-definable text lines
- Input of room and specimen temperatures in °C, °F or K
- Automatic room temperature recording with optional room temperature sensor
- Extensive parameter verification to avoid illegal or inconsistent parameter settings

## • Measurement

- Magnetization and demagnetization curve or demagnetization curve only
- Drift correction: automatic or on command, reminder for drift correction by software if required
- Automatic polarity detection and adjustment of field direction for magnetized rare earth magnets
- Adjustable duration of measurement, usually 20 - 120 seconds depending on specimen type
- Speed enhancement by faster magnetization and sufficiently slow demagnetization
- Measuring speed: constant  $dI/dt$  or  $dJ/dt$  controlled
- Excitation to preset values of field current  $I$ , field strength  $H$  or polarization  $J$
- Automatic stop at  $H_{cJ}$  possible
- Automatic stop at  $-I_{max}$  possible

## • Curves


- Demagnetization curve
- 2- or 4-quadrant hysteresis loop
- Recoil curves or inner (minor) loops, automatically or manual
- $(BH)_{max}$  area
- $BH = \text{const.}$  hyperbolas
- $B = 0$  line

## Evaluation


- Remanence ( $B_r$  or  $J_r$ )
- Coercivity of  $J(H)$  curve  $H_{cJ}$  (intrinsic coercivity)
- Coercivity of  $B(H)$  curve  $H_{cB}$  (normal coercivity)
- Maximum energy product  $(BH)_{max}$
- Maximum field strength  $H_{max}$
- Maximum polarization  $J_{max}$  and  $B_r/J_{max}$  ratio
- $H_k$ : knee shape parameter ( $H$  coordinate of  $J(H)$  curve for  $J = 0.9 \cdot B_r$ )
- $H_x$ :  $H$  coordinate of  $J(H)$  curve for  $J = x \cdot B_r$
- $H_{D5}$  and  $H_{D2}$  results
- Tables of  $J(H)$  and  $B(H)$ , where  $H$  are user defined field strengths
- Tables of  $H(J)$  and  $B(J)$ , where  $J$  are user defined polarizations
- Tables of  $H(B)$  and  $J(B)$ , where  $B$  are user defined flux densities (inductions)
- User values for tables can be predefined and stored. The corresponding functional values are automatically interpolated.
- Output of specimen and measurement parameters and calculated results to ASCII text files, Microsoft Excel® worksheets (xlsx) or Microsoft Excel® xml spreadsheets
- Output of sample and measurement parameters and calculated results to databases
- Temperature correction: conversion of results by means of temperature coefficients of  $B_r$  or  $H_{cJ}$

Microsoft Excel® and a SQL server are not included. Windows 10 normally includes “Microsoft Print to PDF”, which allows an automatic creation of PDFs of the measurement diagrams by *PERMA*, without further user interaction.

## Screenshots

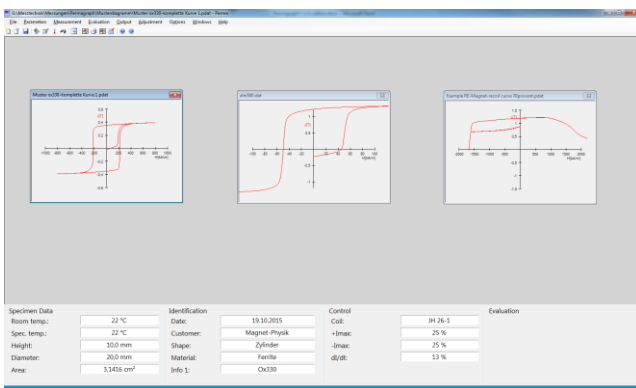


Main window with different measurements

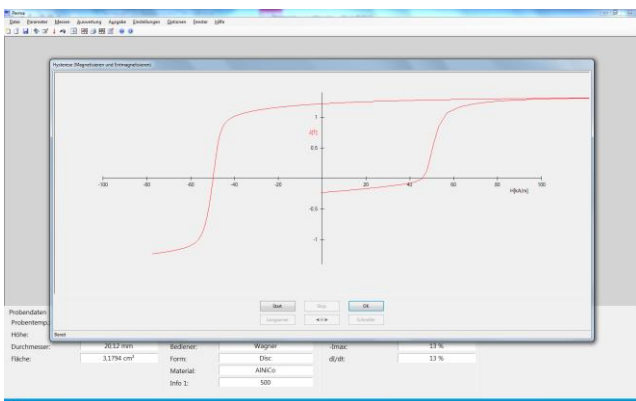

Measurement of a demagnetization curve



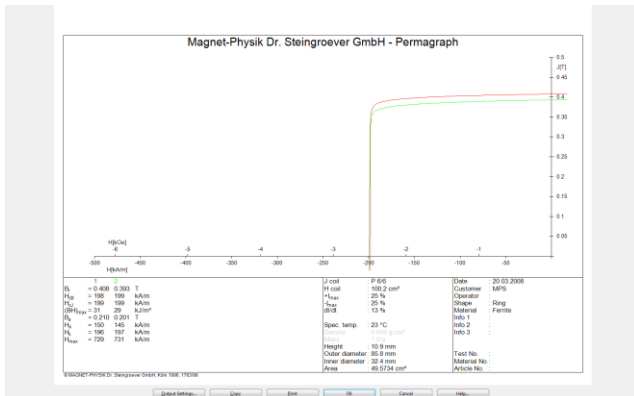

Print preview for multiple curves


## Control

Operation of the PERMAGRAPH® C requires a personal computer to control the measurement course. The computer that is included in the standard packages has the following minimum configuration:

- LCD display
- Keyboard, mouse
- Hard disk or SSD drive
- Network interface
- Interfaces (COM, USB as required)
- Microsoft Windows® operating system (current version)


It is recommended to have the computer configured by MAGNET-PHYSIK. We can only guarantee faultless operation if, apart from the programs installed by us or standard office programs, no further software is installed.

## Example Diagrams

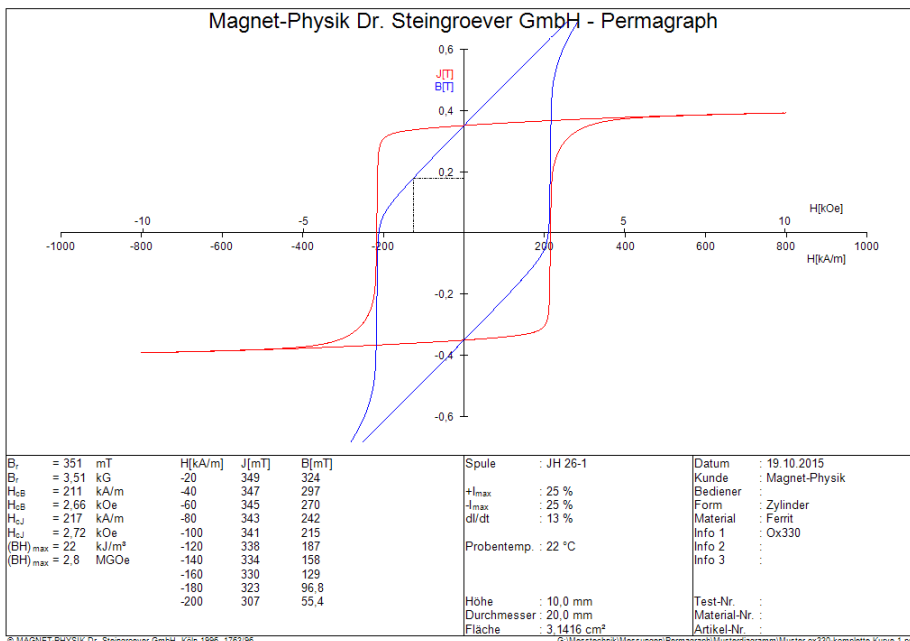


© MAGNET-PHYSIK Dr. Steingroever GmbH, Köln 1986, 1783/96

$J(H)$  and  $B(H)$  hysteresis loops of a ferrite magnet with  $(BH)_{max}$  area. The results and information areas of the diagram are user-configurable.

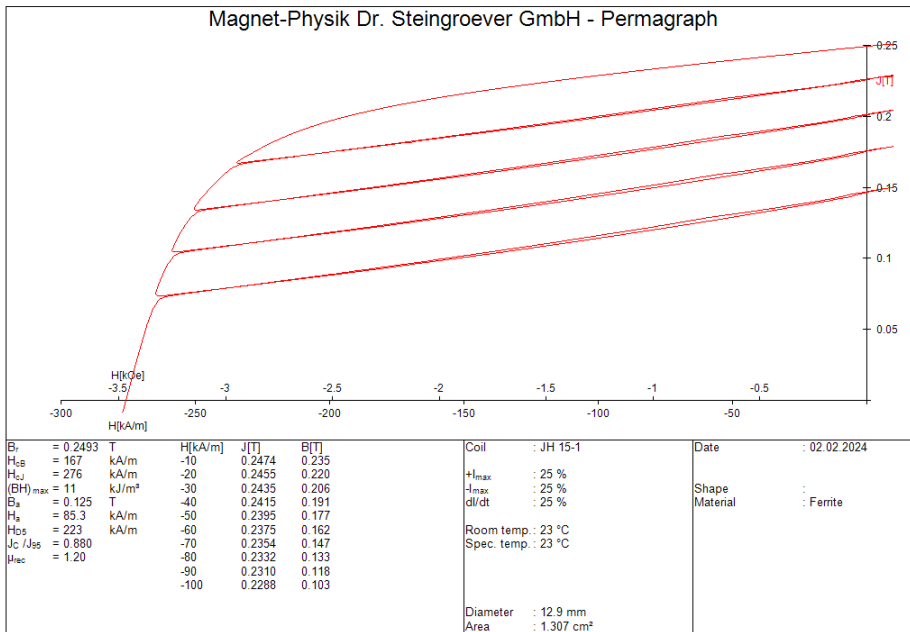


$J(H)$  demagnetization curve of a Ferrite magnet with many recoil loops.

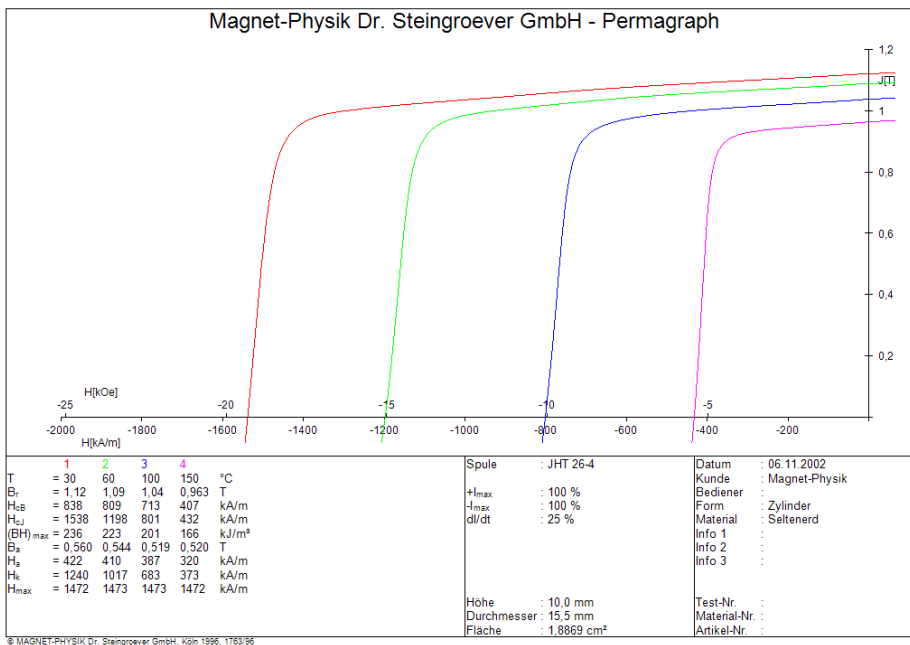

$J(H)$  demagnetization curves of a NdFeB magnet at different temperatures.

## • Electronic Fluxmeter EF 7

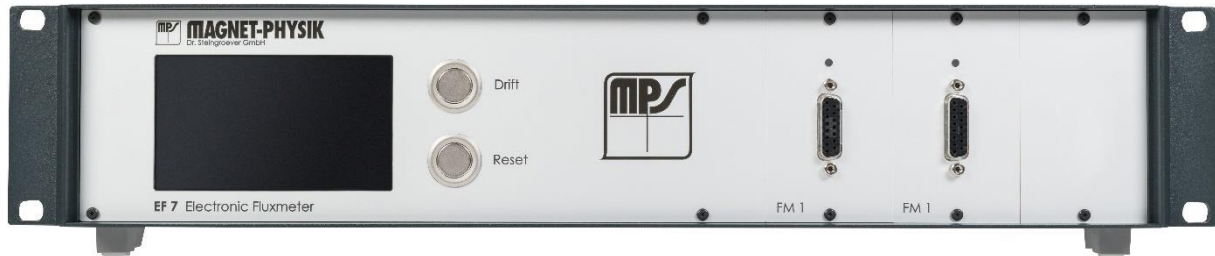


The Electronic Fluxmeter EF 7 is designed to measure the magnetic flux using measuring coils. It comprises a precision digitally compensated DC integrator of high sensitivity and extremely low drift.

For measurements with the PERMAGRAPH® C the fluxmeter is completely controlled from the computer. The measuring coils are equipped with electronic memories containing the coil data. In this way, the connected coils can be automatically identified.

The fluxmeter included with the PERMAGRAPH® C can also be used for other measurement tasks, for example for testing magnets using a moment measuring coil (Helmholtz coil). For details please consult the data sheets "EF 7" and "Measuring Coils for Fluxmeters".

Key features:

- Microprocessor controlled, easy operation
- TFT Touch-Display
- Automatic drift correction
- Unique digitally compensated analog integrator: measuring range limits must not be observed
- Self-calibration by built-in voltage-time reference
- Complete menu control, the most important functions can directly be accessed by function keys
- Automatic calculation of measuring results using the coil parameters
- Directly reading in Volt-Seconds, Weber, Tesla, Gauss or many other units
- Automatic coil recognition and instrument configuration for measuring coils with data memories
- 4 limit comparators with relay outputs for process control
- Modern, compact design
- Directly installable to 19" racks

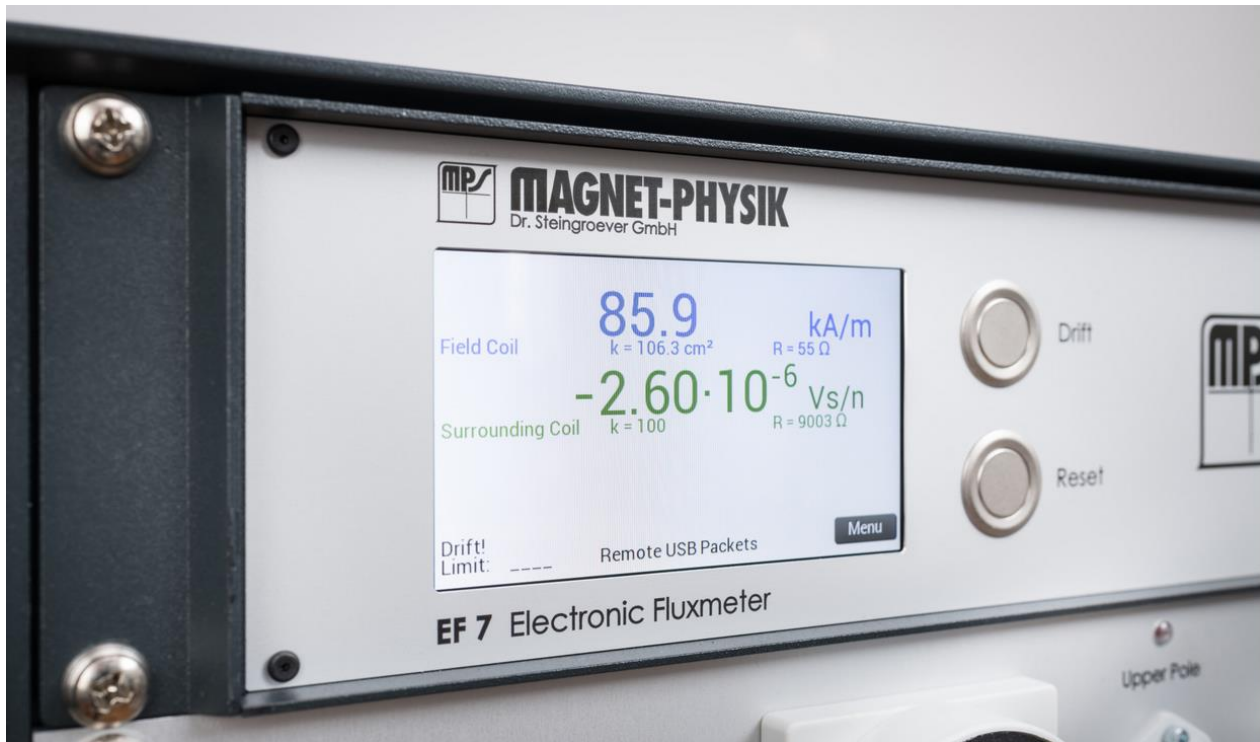

## • Technical Data

Display	TFT 4.3", 95 mm x 54 mm, 480 x 272 points, capacitive touch
Reading	max. 6 digits plus 2 digits for exponent
Resolution	$10^{-4} / 10^{-5} / 10^{-6} / 10^{-7}$ Vs
Upper range limits	none due to digitally compensated integrators
Drift per minute	$< 10^{-6}$ Vs ( $R_i + R_s \geq 10$ k $\Omega$ , $R_s$ = meas. coil resistance)
Basic accuracy	0.25 % of reading
Precision (reproducibility)	0.1 % of reading
Input resistances $R_i$	0 $\Omega$ , 10 k $\Omega$
Maximum input voltage	60 V
Measurements per second	25
Trigger	internal, external
Extreme values	<b> Max.</b> , Max., Min., Max.-Min.
Limit comparator	4 trip points, photo-relay outputs (alternators)

• Electromagnets

### Electromagnet EP 3



High field strength in the air gap: up to 1700 kA/m (21.5 kOe) with 2 mm air gap and straight poles

Exchangeable pole caps with following options:

- Made of soft magnetic steel or FeCo
- Diameter 92 mm straight or tapered
- Without or with pole coils
- Without or with heating (up to 200 °C)
- Air gap plane parallel or with segment profile

Minor heating of the field coils

Easy movement of upper poles by hand wheel

No danger of injuries by skewed yoke frame


Required power supply: 3 kW

### Electromagnet EP 5

The measurement of hysteresis curves of permanent magnets in a measuring yoke requires very high field strengths for the demagnetization of the modern, high coercive materials (Samarium-Cobalt and Neodymium-Iron-Boron).

- Exchangeable pole caps with options as for EP 3
- Easy movement of the upper pole by hand wheel on the front side
- No danger of injuries due to the closed yoke frame
- Required power supply: 3 kW

Achievable flux density in the air gap with iron-cobalt (FeCo) poles tapered to 65 mm diameter for different air gap heights  $h$ :

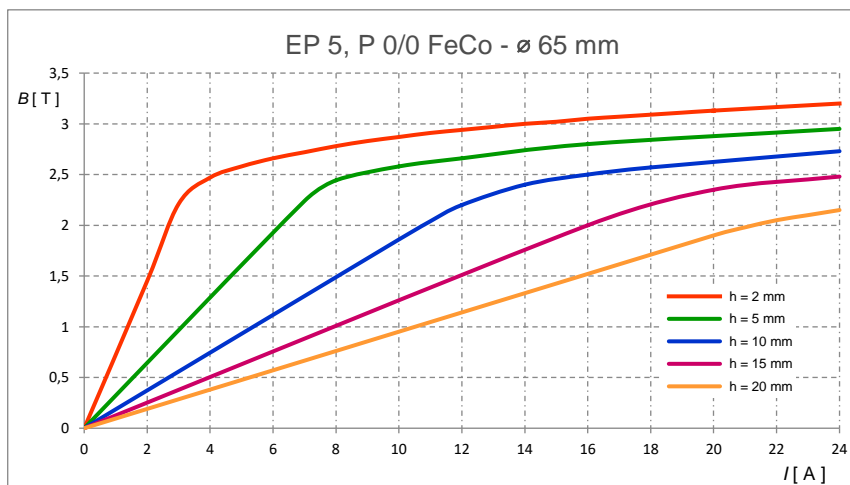

With the electromagnet EP 5, the high field strength in the air gap is achieved by moving the upper field generating coil together with the upper pole in the frame of electromagnet. In this way the magnetic field of the two coils has an optimum effect.

In addition the pole shape and the winding cross sections are of optimum design.

## • Accessories for PERMAGRAPH®

### Electromagnet Poles

Exchangeable pole caps for electromagnets EP 3 and EP 5. Two poles P 0/0 are usually needed. The poles are made of low carbon soft magnetic steel or a (rather expensive) iron cobalt (FeCo, Vanadium Permendur) alloy, which is the material with the highest saturation polarization. Special sizes are available on request.

#### P 0/0 - ø 92 mm

Standard pole for measurement with surrounding coils or flat measuring poles. Made of soft magnetic steel.

#### P 0/0 - ø 65 mm

For measurement with surrounding coils. Tapered from 92 mm to 65 mm. Maximum surrounding coil diameter: 40 mm. Made of soft magnetic steel.

#### P 0/0 FeCo - ø 65 mm

For measurement with surrounding coils. Tapered from 92 mm to 65 mm. Maximum surrounding coil diameter: 40 mm. Made of FeCo.

#### P 0/0 - ø 80 mm

For measurement with surrounding coils or as counterpart to a tapered measuring pole with 80 mm diameter. (Regarding use in an EP 3 please contact MAGNET-PHYSIKER.)



P 0/0 - ø 92 mm



P 0/0 - ø 65 mm  
(EP 5 version)

### Measuring Poles with Pole Coils

With two embedded pole coils for measuring the polarization  $J = B - \mu_0 \cdot H$  at room temperature (suitable for ferrite magnets). Normally pole coils with 6 mm diameter are used. Coils with 3 mm diameter are used to compare the polarization at different locations on the magnet surface.

### Flat Measuring Poles


These measuring poles can be positioned on standard poles P 0/0 - ø 92 mm.

#### P 3/3-15, 92 mm diameter

With two embedded pole coils of 3 mm diameter

#### P 6/6-15, 92 mm diameter

With two embedded pole coils of 6 mm diameter

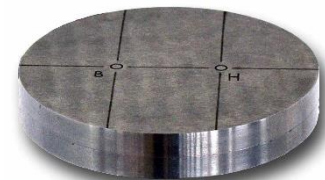

P 3/3-15

### Tapered Measuring Poles

These poles are screwed to the electromagnet. They are tapered from 92 mm to 80 mm. As counterpart a pole P 0/0 - ø 80 mm is required. The poles are made of soft magnetic steel. (Regarding use in an EP 3 please contact MAGNET-PHYSIK.)

#### P 3/3

With two embedded pole coils of 3 mm diameter

#### P 6/6

With two embedded pole coils of 6 mm diameter

#### P 9/9

With two embedded pole coils of 9 mm diameter



P 6/6



• **Accessories for PERMAGRAPH®**

**Field Coil FS 100/2**

For the measurement of the magnetic field strength or flux density on permanent magnets or in magnet systems. Required for the measurement of the field strength  $H$  when measuring poles with pole coils are used.

Thickness: 2 mm, area turns: approx. 100 cm<sup>2</sup>.



**Field Coil FS 100/2**

**Field Coil FS 100/1**

For the measurement of the magnetic field strength or flux density on permanent magnets or in magnet systems. Especially suitable if thin magnet sheets are measured using pole coils.

Thickness: 1 mm, area turns: approx. 100 cm<sup>2</sup>.



**Field Coil FS 100/1**

**J-compensated Surrounding Coils**

With integrated field strength measuring coils for measurements at room temperature. Coil thickness: 1 mm.

**J-compensated Surrounding Coil JH 10-1**

Diameter 10 mm,  
Specimen diameter 5 ... 10 mm

**J-compensated Surrounding Coil JH 15-1**

Diameter 15 mm,  
Specimen diameter 10 ... 15 mm

**J-compensated Surrounding Coil JH 26-1**

Diameter 26 mm,  
Specimen diameter 15 ... 26 mm

**J-compensated Surrounding Coil JH 40-1**


Diameter 40 mm,  
Specimen diameter 26 ... 40 mm

**J-compensated Surrounding Coil JH 60-1**


Diameter 60 mm,  
Specimen diameter 40 ... 60 mm

**Protective covers for JH coils**

For mechanical protection during measurement. The protective cover is clipped onto the coil. This increases the total coil thickness to 7 mm.

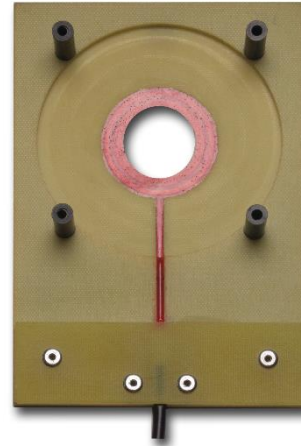

J-compensated Surrounding Coil JH 26-1

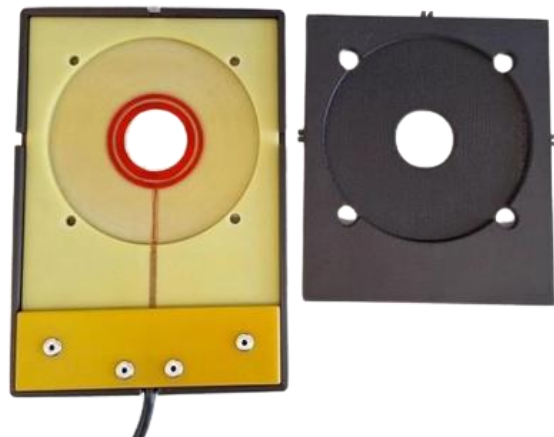

**Other J-compensated surrounding coils, also of rectangular shape, can be supplied on request.**

J-compensated surrounding coils from MAGNET-PHYSIK follow an elaborate design of the windings resulting in separate measuring and compensation coils. This design ensures a real, countable number of turns of the measuring coil. A simpler design approach, which is not used by MAGNET-PHYSIK, requires that an effective number of turns must be determined by measurements on reference specimens, resulting in inferior reproducibility.



## • Accessories for Measurements at Elevated Temperatures

### Heating Poles

For hysteresis measurements at specimen temperatures up to 200 °C. The poles contain a heating element and are tapered from 92 mm diameter to 60 mm diameter. They can be inserted into the electromagnets EP 3 or EP 5. Two poles are needed.

### Heating Poles PT 200

The pole caps are made of low carbon soft magnetic steel.

### Heating Poles PT 200 FeCo

The pole caps are made of a high saturation iron-cobalt alloy. This material provides better measuring accuracy especially in the vicinity of the knee of the demagnetization curve at field strengths above 1000 kA/m.

### Thermocouple TH-J 2

Temperature sensor for the measurement of the specimen temperature, for insertion into heating poles PT 200 or PT 200 FeCo.

### Temperature Surrounding Coils

With integrated field strength measuring coils, for measurements at temperatures up to 200 °C. Coil thickness: 4 mm.

#### Temperature Surrounding Coil JHT 10-4

Diameter 10 mm,  
Specimen diameter 5 ... 10 mm

#### Temperature Surrounding Coil JHT 15-4

Diameter 15 mm,  
Specimen diameter 10 ... 15 mm

#### Temperature Surrounding Coil JHT 26-4

Diameter 26 mm,  
Specimen diameter 15 ... 26 mm

#### Temperature Surrounding Coil JHT 40-4

Diameter 40 mm,  
Specimen diameter 26 ... 40 mm

### Thin Temperature Surrounding Coils

With integrated field strength measuring coils, for measurements at temperatures up to 200 °C. Coil thickness: 2 mm.

#### Temperature Surrounding Coil JHT 10-2

Diameter 10 mm,  
Specimen diameter 5 ... 10 mm

#### Temperature Surrounding Coil JHT 15-2

Diameter 15 mm,  
Specimen diameter 10 ... 15 mm

#### Temperature Surrounding Coil JHT 26-2


Diameter 26 mm,  
Specimen diameter 15 ... 26 mm



Heating Pole PT 200 or PT 200 FeCo



Thermocouple TH-J 2

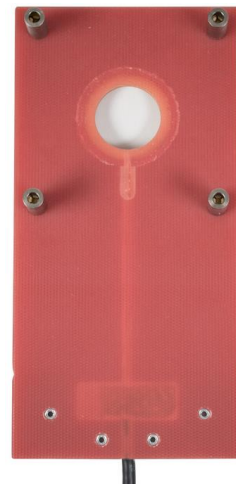

Temperature Surrounding Coil JHT 26-4



Temperature Surrounding Coil JHT 26-2

## • Special Applications

### Segment Poles with Pole Coils

For the measurement of segment magnets. Both the upper and the lower pole contain pole coil. Thus the polarization can be measured either on the lower or the upper side of the segment magnet, or both pairs of pole coils can be connected together to obtain the average.

The set of poles is constructed corresponding to the radii of the segment magnet. Magnets with different sizes and thicknesses but the same radii can be measured with the same pole set.

#### Set of segment poles

Material: soft magnetic steel

#### Set of segment poles

Material: iron-cobalt (FeCo)

The poles are suitable for use in the electromagnets EP 3 and EP 5. For the measurement you also need one field measuring coil FS 100/2 (which is already included in the "Ferrite" package).

Optionally available:

#### Spring Loaded Holding Device TP 2

(not suitable for electromagnets EP 3)

This device makes the mounting of a set of segment poles to the electromagnet faster and easier.

### Segment Poles MC-BLW

For the measurement of flux curves of ferrite segment magnets according to the standard DIN 50472. The set of poles is constructed corresponding to the dimensions of the segment magnet.

A coil in one of the poles senses the total magnetic flux penetrating the magnet. The pole distance is fixed and the measurement is carried out with a well-defined air gap. A sheared hysteresis loop is obtained.

The poles are suitable for the use in the electromagnets EP 3 or EP 5. For the measurement you also need one field coil FS 100/2.



Segment Poles with Pole Coils



Segment Pole MC-BLW

---

## • Combined Instruments

---

### **PERMAGRAPH® - REMAGRAPH® COMBINATION C - 850**

This combination instrument allows the measurement of the static hysteresis loops of soft and hard magnetic materials.

It combines the complete measuring capabilities of PERMAGRAPH® C - 400 and REMAGRAPH® C - 600 into one instrument.

As some components are only required once, the combination offers a reasonable alternative to separate instruments.

Components of the standard package:

- Cabinet for combination C - 850
- Electronic Fluxmeters EF 7-Perma-Rema
- REMAGRAPH® power supply SVR 5
- PERMAGRAPH® power supply SVP 5
- REMAGRAPH® measuring yoke MJR 5 with exchangeable poles
- J-compensated surrounding coil, round JRR
- J-compensated surrounding coil, flat JRF
- Potential coil PS-R-40/58
- Connection box for ring specimens CB-R
- PERMAGRAPH® electromagnet EP 3
- Pole caps P 0/0 92 mm (2 pieces)
- J-compensated surrounding coil JH 26-1 (other diameter on request)
- Computer Hardware
- Control board ST-P/R 2
- Software PERMA and REMA
- Introduction and training in our premises

All accessories of the PERMAGRAPH® C can also be used with the combination.

You can find a detailed description of all features of the REMAGRAPH® C and of the optional accessories in the individual specification of this instrument.

## • Further Accessories

### Nickel Samples

Cylinders with a height of 10 mm and cross-sectional areas of 0.5 cm<sup>2</sup>, 1.0 cm<sup>2</sup>, 2.0 cm<sup>2</sup> or 5.0 cm<sup>2</sup> are available. A proprietary calibration certificate is included. Saturation polarization: 0.6 T ± 2 %.

Nickel samples are traditional references for calibrating the measurement of *J* but they are not necessary for operation of the PERMAGRAPH®.



Nickel Sample

### Reference Magnet

Isotropic ferrite magnet embedded in a brass frame, diameter 24 mm, for testing and calibrating the PERMAGRAPH®, with proprietary calibration certificate and measuring diagram.



Reference Magnet

### Room Temperature Sensor TS-USB

Temperature sensor for USB connection. Using this sensor the room temperature can be shown on the computer display and automatically be saved with the measurement.



TS-USB

## • Services

### Taking into Operation and Training – at the Premises of MAGNET-PHYSIK

Training in the operation and software of the computer controlled PERMAGRAPH®.

The training takes 1 day and is included in the standard packages. The customer bears all costs in connection with the journey, stay, accommodation, etc.

### Taking into Operation and Training – at the Premises of the Customer

Training in the operation and software of the computer controlled PERMAGRAPH®.

The training takes 1 day and is at the customer's expenses. Additionally, the customer bears the costs for our employee in connection with the journey (incl. expenses for travel time, stay, accommodation, etc.).

*Due to continuous product improvements, specifications are subject to change without notice.*

### MAGNET-PHYSIK Dr. Steingroever GmbH

Emil-Hoffmann-Straße 3, 50996 Cologne, Germany  
Phone: +49 2236 3919-0 • Fax: +49 2236 3919-19  
info@magnet-physik.de  
www.magnet-physik.de

### MAGNET-PHYSICS Inc.

6330 East 75th Street, Suite 224, Indianapolis, IN 46250, USA  
Phone: +1 317 577 8700 • Fax: +1 317 578 2510  
info@magnet-physics.com  
www.magnet-physics.com